


LightScanner® Master Mix

Information Sheet

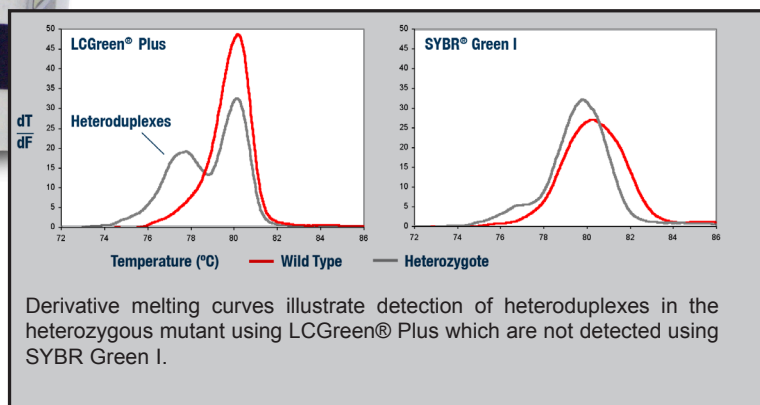

Designed for Hi-Res Melting® Applications

LightScanner Master Mix brilliantly provides a pre-optimized solution specifically designed to deliver superior PCR performance for Hi-Res Melting applications on the LightScanner Instrument.

Product Benefits


- **Reduce time, save money**
easy-to-use premixed formula
- **Secure downstream success**
go directly from PCR to melting curve analysis, eliminating the risk of contamination.

Idaho Technology is now



Superior PCR Performance

The two images below show the amplification of the Hepatic Lipase gene (301 bp) using 100 ng human genomic DNA using the LightScanner Master Mix. Superior reproducibility is demonstrated by the overlapping curves of 24 replicate samples of both the wild type and mutant samples.

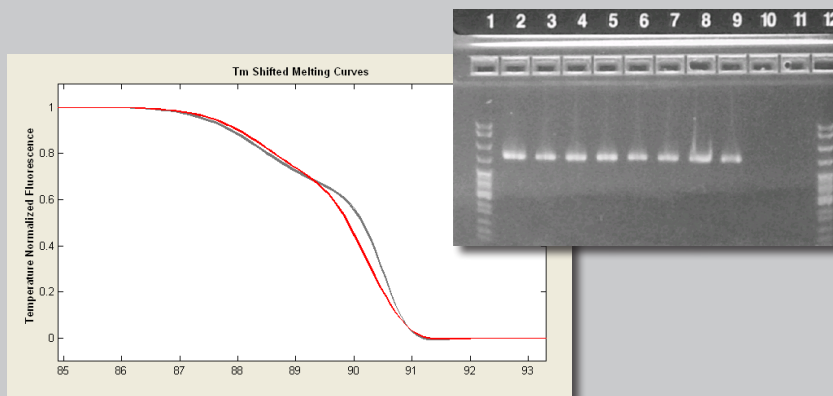


LightScanner® Work Flow

The LightScanner System was developed to integrate seamlessly into existing laboratory work flows. The process is nondestructive so samples can be recovered for additional analysis.


Step 1: Wet Reaction Setup

Follow basic PCR reaction setup using the Master Mix, primers and DNA. The Master Mix contains LCGreen® Plus.

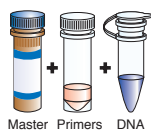


Step 2: Place Mix in Plate with Oil Overlay

Preload PCR plate with mineral oil, add wet mix, seal with film and centrifuge briefly at 2500 rpm.

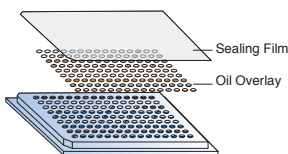

Step 3: Thermocycle Samples

Follow the recommended protocols to perform PCR.

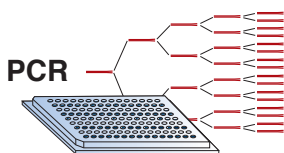


Step 4: Melt Samples in LightScanner

Following PCR, insert the 96- or 384-well plate into a LightScanner to melt the samples.




Step 5: Analyze Melt Data

Following the melt, use the LightScanner software to manage and analyze the data.

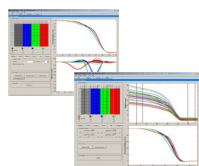

Step 6: Scanning is Nondestructive

Samples can be recovered for additional analysis, sequencing, gel electrophoresis, remelting, etc.

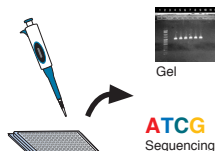

Approximately
5 to 8 Minutes

Shipping and Storage

- Store at -20 °C for up to 1 year
- Once thawed store at 4 °C for 2 weeks
- LightScanner Master Mix is shipped on dry ice

Kit Contents

| Quantity | Description |
|-------------|--------------------------------------|
| 1 x 100 Rxn | 2.5 X Master Mix |
| 1.5 mL | 10 mM MgCl ₂ |
| 1.5 mL | Reagent Grade Water |
| 1 | LightScanner Master Mix User's Guide |

Ordering Information

| No. of Rxns | 100 | 500 | Larger sizes Inquire |
|-------------|---------------|---------------|----------------------|
| Catalog No. | HRLS-ASY-0002 | HRLS-ASY-0003 | |

BIOKÉ
sharing knowledge

Plesmanlaan 1d
2333 BZ Leiden
The Netherlands
T. +31 (0)71 568 10 00
T. Belgium: 0800 71640
F. +31 (0)71 568 10 10
info@bioke.com
www.bioke.com

The purchase of the LightScanner Master Mix includes a limited, nontransferable license, under specific claims of one or more U.S. patents as listed on BioFire Diagnostics, Inc.'s Web site (<http://www.biofiredx.com/LegalNotices/>) and owned by the University of Utah Research Foundation and/or BioFire Diagnostics, Inc., to use only the enclosed amount of product according to the specified protocols. No right is conveyed, expressly, by implication, or by estoppel, to use any instrument or system under any claim of such U.S. patent(s), other than for the amount of product contained herein. Purchase of the LightScanner Master Mix does not convey any PCR license. The LightScanner Master Mix is licensed under U.S. Patent Nos. 5,338,671 and 5,587,287. The purchase of the LightScanner Master Mix includes a limited, non-transferable license for all fields other than human or veterinary in vitro diagnostics under U.S. Patent No. 5,871,908, owned by Evo-tec Biosystems GmbH and licensed to Roche Diagnostics GmbH, to use only the enclosed amount of product according to the specified protocols. No right is conveyed, expressly, by implication, or by estoppel, to use any instrument or system under any claim of U.S. Patent No. 5,871,908, other than for the amount of product contained herein.

LCGreen, Hi-Res Melting, LightScanner are trademarks of BioFire Diagnostics, Inc. LightCycler is a trademark of a member of the Roche Group. SYBR is a trademark of Molecular Probes.

References

1. Olsen RK, Dobrowolski SF, Kjeldsen M, Hougaard D, Simonsen H, Gregersen N, Andresen BS. (2010). **High-resolution melting analysis, a simple and effective method for reliable mutation scanning and frequency studies in the ACADVL gene.** J Inher Metab Dis. 33(3): 247-60.
2. Hulvey J, Young J, Finley L, Lamour K. (2010). **Loss of heterozygosity in Phytophthora capsici after N-ethyl-nitrosourea mutagenesis.** Mycologia. 102(1): 27-32.
3. Audrezet MP, Dabricot A, Le Marechal C, Ferec C. (2008). **Validation of high-resolution DNA melting analysis for mutation scanning of the cystic fibrosis transmembrane conductance regulator (CFTR) gene.** J Mol Diagn. 10(5): 424-34.
4. Wittwer CT, Reed GH, Gundry CN, Vandersteen JG, Pryor RJ. **High-resolution genotyping by amplicon melting analysis using LCGreen.** Clin Chem. 2003 Jun; 49(6):853-60.

Technology Comparisons

1. Chou, L.-S., Lyon, E. and Wittwer, C. T. **A comparison of high-resolution melting analysis to denaturing high performance liquid chromatography for mutation scanning: cystic fibrosis transmembrane conductance regulator gene as a model.** Am J Clin Pathol, 2005 Sep; 124(3):330-8.
2. Herrmann MG, Durtschi JD, Bromley LK, Wittwer CT, Voelkerding KV. **Amplicon DNA melting analysis for mutation scanning and genotyping: Cross-platform comparison of instruments and dyes.** Clin Chem. 2006 Mar; 52 (3):494-503. Epub 2006 Jan 19.