



NextGENe[®]

Network License Renewal Guide

1. Open the SoftGenetics License Server Manager that is running on Server Computer. You can open this service by right-clicking on the  icon from icon tool tray and selecting "Show".
2. To view licensing information, click on "Licensing Information" from the View menu. When Valid Period is less than 1 day, the license is expired; license renewal and registration is required.


#	Product Name	Max Number	Current Number
1	NextGENe	2	0
2	NextGENe Viewer	8 (additional)	0
3	Mutation Surveyor	2	0
4	GeneMarker HID	5	0
5	ChimerMarker	5	0

You have registered:

Product Name: NextGENe
Valid period: 0 days
Max user: 2
Current user: 0

Close

1. Register NextGENe by clicking on “Register” from Help menu.
2. Select “NextGENe” as the Product.
Input your Account, Password and Email and click Register.
 1. Upon successful registration, you will be prompted to restart License Server Manager for the changes to take effect.
 2. If unsuccessful, click on the Offline Registration tab.

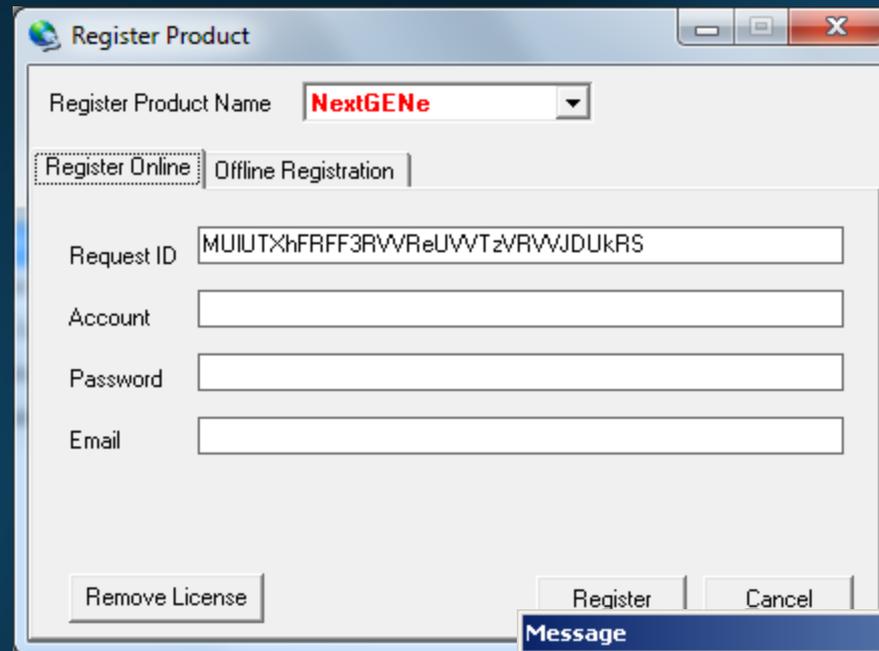

Register Product

Register Product Name **NextGENe**

Register Online | Offline Registration

Request ID MUIUTXhFRFF3RVVReUWWTzVRWJDUKRS

Account

Password

Email



Remove License Register Cancel

Message

The SoftGenetics License Server Manager needs to be restarted for changes to take affect. Please make sure that no clients are connected before restarting.

Restart Now Restart Later

1. Email the request ID and your account credentials to tech_support@softgenetics.com, and you will received your Registration ID.
2. Paste Registration ID into the Register ID field and click Register.
3. You will be prompted to restart License Server Manager for the changes to take effect.

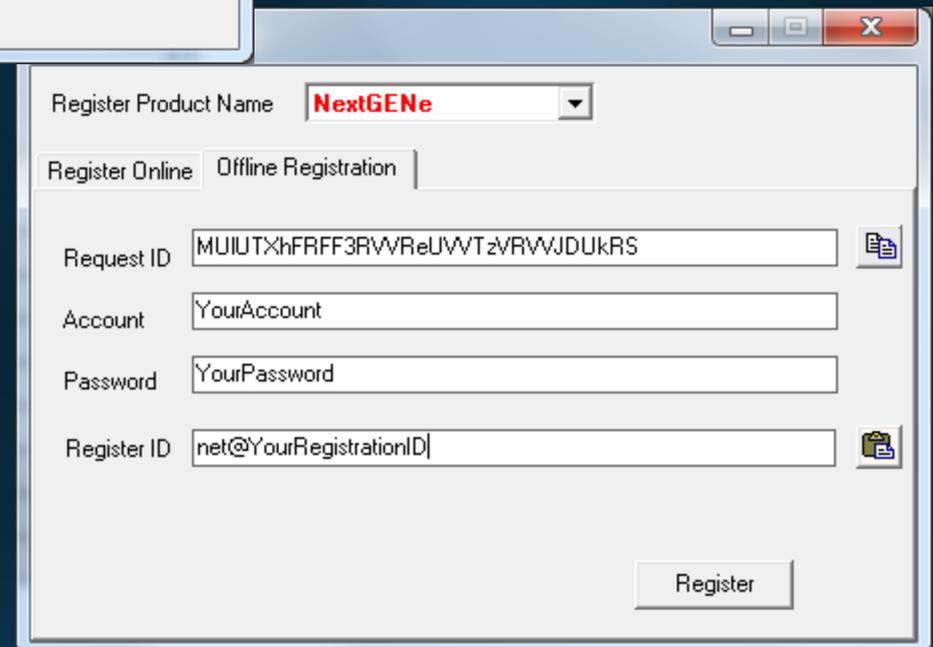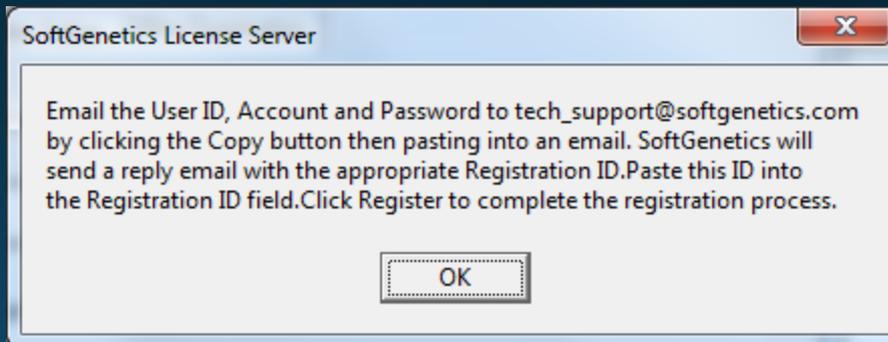

NextGENe will now open with full functionality from all client computers.

