

**TUTORIAL FOR THE CORRECT USE
OF DISPOSABLE HOMOGENIZATION CONSUMABLES**

Description:

disposable devices for homogenization of biological samples for molecular extractions

Maximum volume
(sample + solvent)

■ **NEVER USE CONSUMABLES DRY**

■ **SEAL CONSUMABLES COVER BEFORE PUTTING INTO MEDIMACHINE II**

Fragmenting the sample as much as possible

Avoid too heavy or compact samples that block the start of the rotation of the blades


Sample withdrawal:

FOODcon7: completely remove the cover and withdraw the liquid homogenate

Prypcons: open the transparent cover by sliding it and withdraw the liquid homogenate

FOODcon7

Prypcons

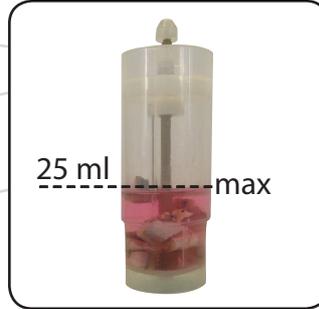

FOODcon7
&
Prypcons