



Instructions for Use

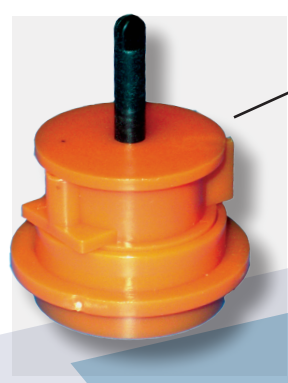


Cod. 79300

Ideal volume
1 ml

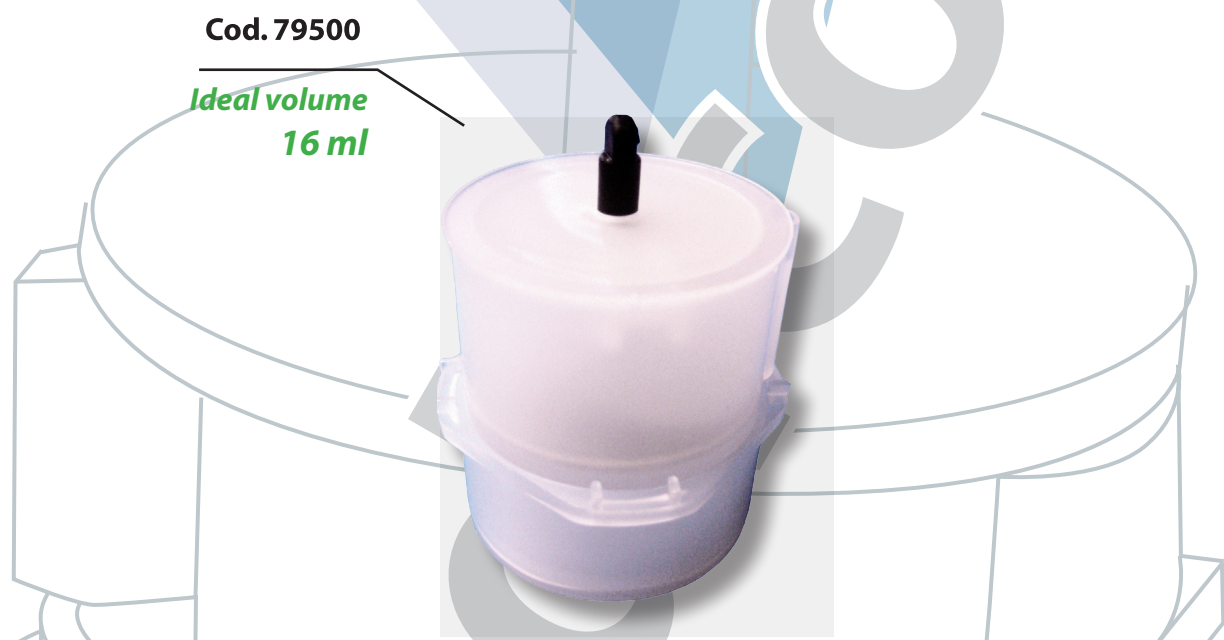

Cod. 79500

Ideal volume
16 ml

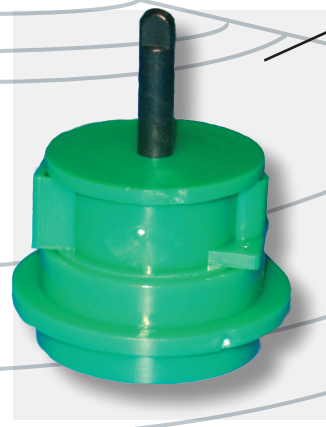

Cod. 79400

Ideal volume
1 ml




www.ctsv.biz



Description:


Disposable devices for the dissociation of biological samples for cells extraction

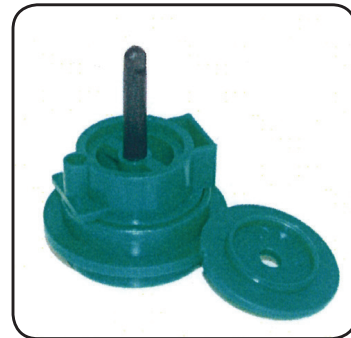

Too much liquid please check is the sample floating.
(The sample is not grinded)

Maximum volume of liquid to be added to the sample = 1 ml


Not enough liquid will leave the grater dry
and can possibly damage the sample

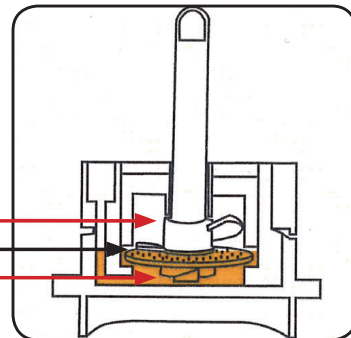

Ideal volume Medicons 79500 = 16 ml

Cut the sample into small pieces and place it on grater

Avoid placing the sample on the motor.

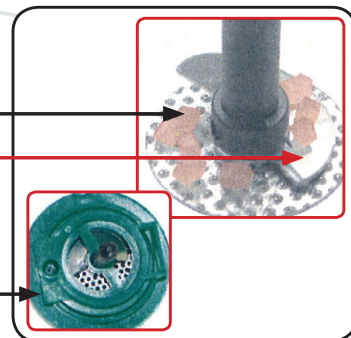

Lift the lid before inserting the Medicons into the Medimachine
and start operation as instructed.


SAMPLE RECOVERY

The liquid containing the cell suspension is collected
through the opening using a

NEEDLELESS SYRINGE



cell suspension

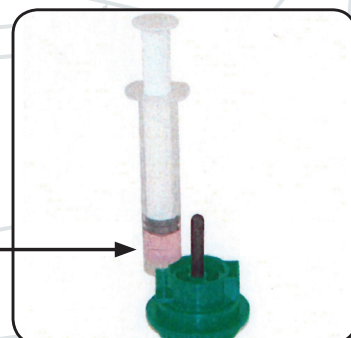

DO NOT DISASSEMBLE MEDICONS

Procedure valid for all types of Medicons